


# General Title: VPE Software: SecureVPE Suggested Access Level Table

<b>Abstract:</b>	This document provides details of suggested security parameters inside the SecureVPE application.
<b>Applicability:</b>	This article applies to the SecureVPE and QuickVCA applications which are a part of the C Technologies, Inc. SoloVPE Software.
<b>Symptom:</b>	N/A
<b>Cause:</b>	Use of the SecureVPE Software and the company's need to configure the software to best meet their internal requirements. It is the responsibility of each company to implement this software and environmental requirements relating to file/data integrity.

## Detailed Info: Variable Pathlength Security Guidance for Windows NTFS Structure

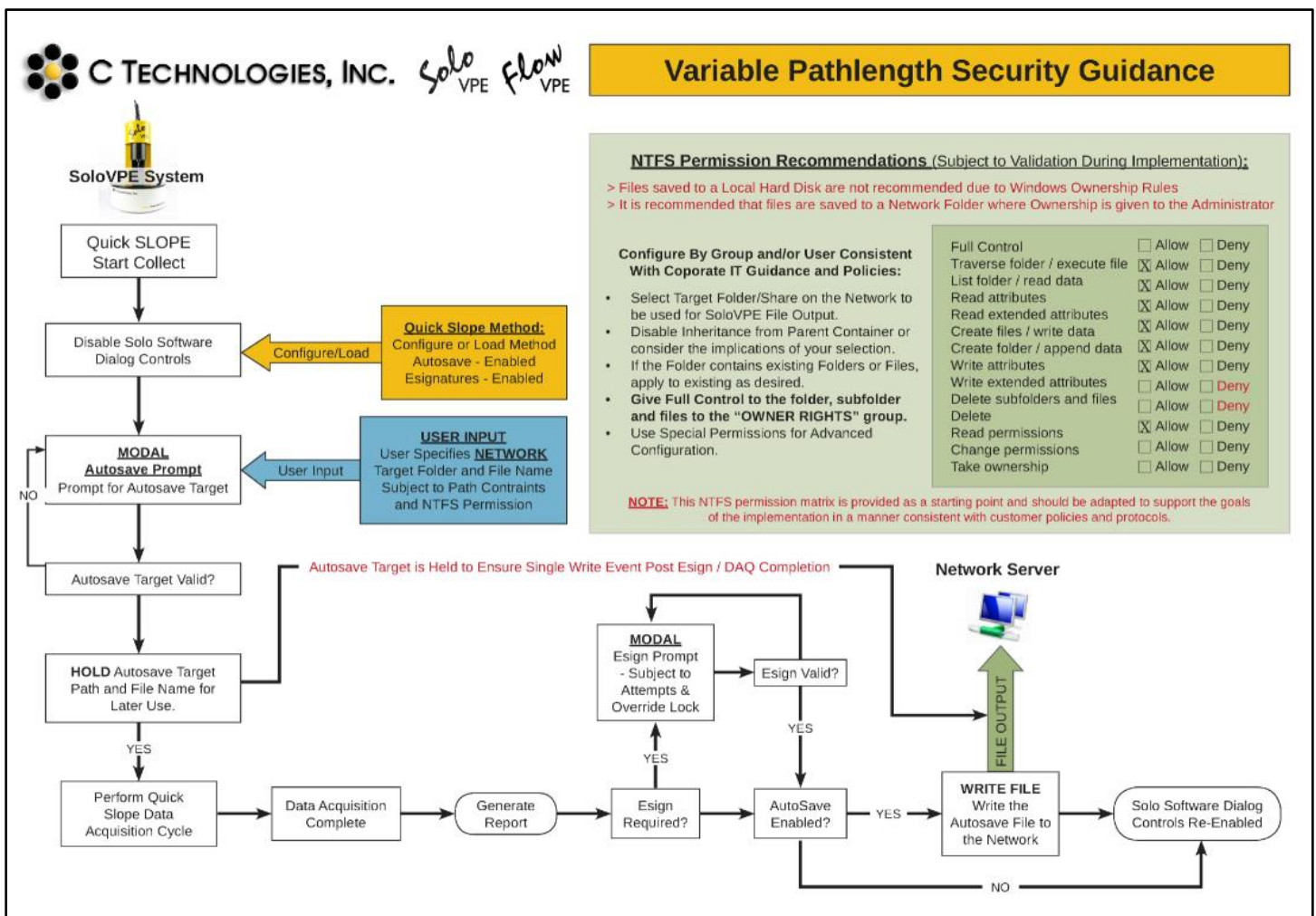

## Detailed Info: Variable Pathlength Security Guidance SecureVPE for SoloVPE Software

<u>Group</u>	<u>Privilege</u>	<u>Description</u>	VPE Admin	VPE Power User	VPE User	VPE Reviewer
<b>SoloVPE</b>						
<b>Application</b>	SoloVPE Software Access	Authority to Run/Use Application	x	x	x	x
	SoloVPE Administration Access	Authority to Run/Use Application	x			
	SecureVPE Access	Authority to Run/Use Application	x			
<b>Method</b>	Method Modification Rights Enabled	Authority to Modify Methods	x	x		
	Advanced Setting Access Enabled	Authority to Access Advance Setting	x	x		
	DAQ View Enabled	Authority to Access DAQ View (open method only)			x	
	DAQ View Optional	Authority to Access DAQ View (optional)	x	x		
	Quick Methods Enabled	Authority to Add Quick Method	x	x		
	Force Disable Quick Slope Alerts	Authority to Disable Quick Slope Alerts				
<b>Sidebar</b>	Report Sidebar Button	Report Sidebar Button Visible	x	x		
	Manual Sidebar Button	Manual Sidebar Button Visible	x	x		
	Section Sidebar Button	Section Sidebar Button Visible	x	x		
	Analyze Sidebar Button	Analyze Sidebar Button Visible	x	x		
	Quick Survey Sidebar Button	Quick Survey Sidebar Button Visible	x	x		
	Quick Slope Sidebar Button	Quick Slope Sidebar Button Visible	x	x	x	
	EC Library Sidebar Button	Ext Coefficient Library Sidebar Button Visible	x	x		x
	Home Sidebar Button	Home Sidebar Button Visible	x	x	x	
	New Run Sidebar Button	New Run Sidebar Button Visible	x	x	x	
	Save Data Sidebar Button	Save Data Sidebar Button Visible	x	x	x	x
<b>Toolbar</b>	Factor Scale Toolbar Button	Factor Scale Toolbar Button Visible	x	x		
	Clear Rectangles Toolbar Button	Clear Rectangles Toolbar Button Visible	x	x		
	Slope Tool Toolbar Button	Slope Tool Toolbar Button Visible	x	x		
	Arrange Graphs: HB	Arrange Graph HB Visible	x	x	x	
	Arrange Graphs: HT	Arrange Graph HT Visible	x	x	x	
	Arrange Graphs: VR	Arrange Graph VR Visible	x	x	x	
	Arrange Graphs: VL	Arrange Graph VL Visible	x	x	x	
	Volume Calculator Toolbar Button	Volume Calculator Button Toolbar Visible	x	x		
	Manual eSignature Toolbar Button	Manual eSignature Toolbar Button Visible	x	x	x	
	Quick Check Toolbar Button	Quick Check Toolbar Button Visible	x	x	x	
	Plugins Enabled	Additions or Patches to the SoloVPE Software (inactive)				
<b>Report</b>	Report Modification Rights Enabled	Authority to Modify All Reports Enabled				
	Quick Slope Report Options Enabled	Ability to use Quick Slope Report Options (Advanced Tab)	x	x		
<b>Administration</b>	Measurement Continuation Allowed	Continue Performing Sample Measurement	x	x	x	
	Set Zero Allowed	Ability to Set Zero Position	x			
	Quick Check Administrator Access	Authority to Access Quick Check Administrator	x			
	Optimize Coupler Access	Authority to Access Optimize Coupler	x			
	Set Zero Wizard Access	Authority to Access Set Zero Wizard	x			
	Manual Limit Override Access	Authority to Access Manual Limit Override	x			
	EC Edits Allowed	Authority to Allow Edits in EC Library	x	x		
	Allow Default Path Subfolders	Authority to Save Files in Subfolders within the Applicable Default Folder	x*	x*	x*	x*
	Force Data Save Default Path	Controls Save Data Events to the Default Paths Configured in the SoloVPE Admin Software.	x	x	x	x
	Force Method Save Default Path	Controls Save Method Events to the Default Pathing Configured in the SoloVPE Admin Software.	x	x	x	
	Coupler Check Access	Authority to ccess the Coupler Check in the Quick Check Application	x	x	x	
<b>Autosave</b>	Autosave Optional	Authority to Autosave Optional				
	Incremental Autosave Enabled	Authority to Incremental Autosave Enabled	x	x	x	
	Force Autosave Default Path	Authority to Force Autosave Default Path	x	x	x	

	Prevent Autosave to Local Hard Drive	Authority to Prevent Autosave to Local Hard Drive	x	x	x	
<b>ESign</b>	ESign Override Rights Enabled	ESign Override Rights Enabled	x			
	ESignatures Enabled	ESignatures Enabled	x	x	x	x
	ESign Quick Check Event Optional	Authority to eSign Quick Check Event	x	x	x	
	Post Manual eSign Save Required	Post Manual eSign Save Required if Checked	x	x	x	
	ESign Pre-Save Event Optional	Authority to eSign Pre-Save Event	x	x	x	
	ESign Quick Methods Optional	Authority to eSign Quick Methods	x	x	x	
	ESign EC Events Optional	Authority to eSign EC Events	x	x		
<b>Security</b>	Inactivity Timer Enabled	Inactivity Timer Enabled (while in software)	x	x	x	x
	Perimeter Authentication Enabled	Perimeter Authentication Enabled (prior to software)	x	x	x	x
	Admin Override Rights Enabled	Authority to Override Administrative Rights	x			

\*It is recommended that all folders only be created by IT/network administrators. Users should not save data to folders created by non-administrators to avoid ownership privileges.

### Detailed Info: Variable Pathlength Security Guidance SecureVPE for QuickVCA Software

<u>Group</u>	<u>Privilege</u>	<u>Description</u>	VPE Admin	VPE Power User	VPE User	VPE Reviewer
<b>QuickVCA</b>						
<b>Application</b>	QVCA Software Access	Authority to Run/Use application	x	x	x	x
<b>ESign</b>	QVCA eSign Enabled	ESignatures Enabled for QVCA Software	x	x	x	x
	QVCA Post Manual eSign Save Required	ESignatures Enabled for QVCA Software	x	x	x	
<b>Method</b>	QVCA Method Modification	Authority to Modify Methods in QVCA	x	x		
<b>Report</b>	QVCA Report Modification Rights Enabled	Authority to Report Modification				
<b>Autosave</b>	QVCA Autosave Optional	Authority to Autosave Optional				
	QVCA Incremental Autosave Enabled	Authority to Incremental Autosave Enabled	x	x	x	
	QVCA Force Data Save Default Path	Authority to Force AutoSave Default Path	x	x	x	
	QVCA Prevent Autosave to Local Hard Drive	Authority to Prevent Autosave to Local Hard Drive	x	x	x	x
<b>Security</b>	QVCA Perimeter Authentication Enabled	Perimeter Authentication Enabled	x	x	x	x
<b>Administration</b>	QVCA Allow Default Path Subfolders	Authority to Force Default Path Subfolders	x	x	x	x
	QVCA Force Data Save Default Path	Authority to Force Save Default Path	x	x	x	x
	QVCA Force Method Save Default Path	Authority to Force Method Save Default Path	x	x	x	


**Document Info: KB17007**

<b>Revision History</b>			
<b>Rev</b>	<b>Date</b>	<b>Changes</b>	<b>Initials</b>
00	2018-02-28	Initial Release	JF
01	2018-12-14	Addition of SecureVPE Check Points	SH
02	2019-07-16	Addition of SecureVPE Check Points/Addition of VPE Reviewer	NH

**Prepared By:**

**C Technologies, Inc.**

685 US Route 202 Bridgewater, NJ 08807

 (908) 707-1009

 [support@ctechnologiesinc.com](mailto:support@ctechnologiesinc.com)