


Packaging Instructions for OPUS® 2.5 - 8 Pre-packed Columns

1. Remove OPUS® Pre-packed Column(s) from storage location.
2. Gather OPUS® Pre-packed Column shipping assembly components ([Figure 1](#)):
 - Individual product box (white)
 - Foam inserts (x2)
3. Assemble individual OPUS® Pre-packed Column box (white; [Figure 1](#)) and set aside.
4. Secure one foam insert onto each end of OPUS® Pre-packed Column.
5. Place OPUS® Pre-packed Column into individual product box:
 - a. Position individual product box such that the longest face is parallel to the unpackaging table/bench.
 - b. Unfold top flap of individual product box.
 - c. Slide foam inserts and OPUS® Pre-packed Column into box as one cohesive unit.
6. Fold down top flap down to secure OPUS® Pre-packed Column within individual column box.
7. If shipping multiple OPUS® Pre-packed Columns with 2.5 - 8 cm ID, repeat Steps 1 - 6 for each column.
8. Place each individual product box within a larger outer box for shipping.
9. Affix address label and packing list to one side of the box.
 - If applicable, affix Refrigerate Upon Arrival stickers.
10. For international shipments, attach the Commercial Invoice to the outside of the box
11. Bring box/es to shipping pick-up area.
12. If OPUS® Pre-packed Column is being shipped to Repligen, contact Complaints@repligen.com with the following information:
 - Part number
 - Lot number
 - Serial number
 - Associated PO number
 - Associated RMA number (provided by Repligen Customer Service)
 - Customer contact name
 - Weight and dimensions of final packaging
 - Image of final packaging
 - MUST include picture to trigger pickup of shipment

Figure 1. OPUS® Pre-packed Column individual product box, 2.5 - 8 cm ID

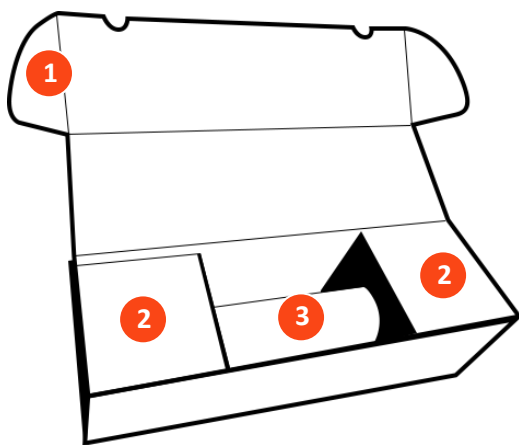

1. Individual product box (white)
2. Foam insert (x2)
3. OPUS® Pre-packed Column

OPUS® Pre-packed Chromatography Columns User Guide



Full User Guide available at
www.repligen.com/resources/quality

Contact information

For more information or questions, please contact Repligen:

Customer Support USA

customerserviceUS@repligen.com

+1-781-250-0111

Repligen Corporation

41 Seyon Street Building #1, Suite 100
Waltham, MA 02453

Customer Support Europe

customerserviceEU@repligen.com

+31 76 5719 419

Repligen Europe B.V.

Hoeksteen 13
4815 PR BREDA
The Netherlands

©2021 Repligen Corporation. All rights reserved. The trademarks mentioned herein are the property of Repligen Corporation and/or its affiliate(s) or their respective owners.